




TOX For Chlorinated Sites

Please read all the instructions before sampling to ensure proper test results.

Additional TOX Sampling Steps

1. Collect sample in bottle so that it is nearly full (to bottom of the neck of the container). **Do not overflow!**
 - We advise you to wear safety glasses and gloves.
2. Add entire contents of small clear vial labeled Sulfuric Acid to the sample bottle. **Do not overflow!**
 - Caution: Avoid skin contact! If acid is spilled or comes in contact with skin, wash thoroughly with water.
3. Fill bottle with sample so that no air remains when capped.
4. Indicate sampling date, time, site and name of sampler on the bottle and on the enclosed Chain of Custody form. Information on the Chain of Custody and labels must match and be complete.

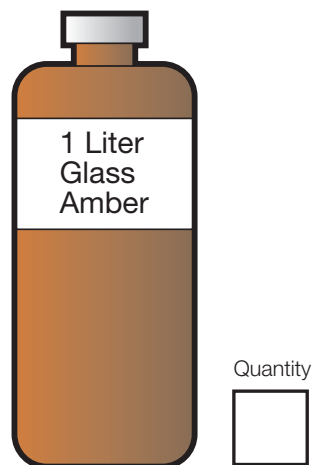

Call 1.574.233.4777 or 1.800.332.4345 or visit us at www.ul.com/water
UL, 110 S. Hill Street, South Bend, IN 46617